**Additional information**  
**Zusatzinformation**  
**Informations complémentaires**  
**Información adicional**

**Additional information**  
**Model PG43SA-S with vent valve**

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Prior to starting any work, read the operating instructions!  
Keep for later use!

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### Supplementary documentation:

- ▶ This additional information applies in conjunction with the operating instructions “Pressure gauge for sanitary applications, models PG43SA-S, PG43SA-C” (article number 11532319).

**This additional information supplements the respective chapters of the above-mentioned document.**

## 2. Design and function

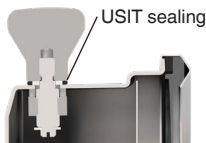
### 2.3 Vent valve

By opening or closing the vent valve, the case will be vented to or sealed from the atmosphere.


Vent valve

Vent valve closed

Vent valve open

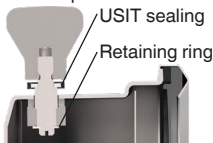- To close: Turn approx. 3 turns in the direction of the “CLOSE” arrow, until the USIT sealing is pressed together fully.
- To open: Turn approx. 3 turns in the direction of the “OPEN” arrow, until the retaining ring prevents the vent valve from being unscrewed further and thus lost.

For any external cleaning (“wash-down”), e.g. in a cabinet washer, the vent valve must be closed.

## 5. Commissioning, operation / 9. Specifications

### 5. Commissioning, operation

EN

#### 5.6 Commissioning and operation

Open the vent valve for pressure compensation and leave it open during the measurement in order to ensure a correct pressure display. The model PG43SA-S with vent valve has no vent hole.

#### 5.7 External zero point setting

External zero point setting is not available for the model PG43SA-S with vent valve.

#### 5.8 Autoclaving

The following is valid only for the model PG43SA-S with the option “Autoclavable version”, identifiable through the dial marking “AUTOCLAVABLE”.

To carry out the autoclaving process correctly, the vent valve must be open.

### 7. Maintenance and cleaning

#### 7.2 Cleaning

For any external cleaning (“wash-down”), e.g. in a cabinet washer, the vent valve must be closed. This is needed in order to prevent the cleaning agent finding its way into the case interior.

### 9. Specifications

#### 9.6 Ingress protection per IEC/EN 60529

- IP68 (with closed vent valve)
- IP54 (with open vent valve)

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All rights reserved. / Alle Rechte vorbehalten.
WIKA® is a registered trademark in various countries.